Connecting your FH III to your PC to use Hunter III Software


Please read through before attempting so you know what to do

1. Plug USB – RS232 converter (only converter not connected to the FH III) into your PC



2. Go to Device Manager on your PC


A. Right click on properties Match comm settings in the FH III (under communications setup).

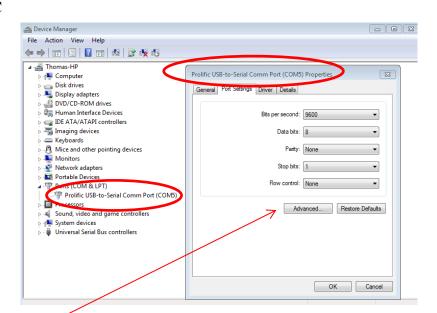

B. If you have issues click on advance – (Under properties)

Now lower the TX and RX buffer to the lowest number to correct for communication errors.

Click ok to save... May have to hit OK again to save under properties also.

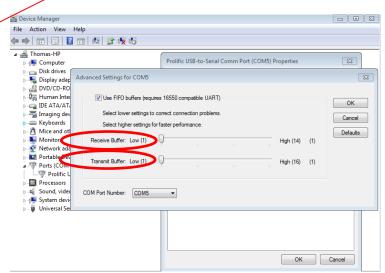


3. Match comm settings in Device manager and the FH III (under communications setup)






- 4. Turn off the FH III, wire up the RS232 wires.
- A. Black to TX
- B. Red to RX

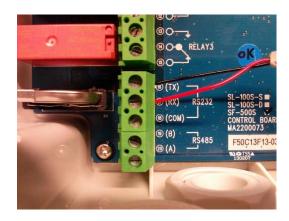

- 5. Now power on your FH III, plug the USB adaptor to the RS232 wire connector.
- 6. Click on

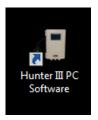

7. Go to File then click on Acquire.





8. Check your comm port and baud rate. (make sure they match your previous comm Settings).

Click Acquire... And now wait for the download.

